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```
$ sh ./topspin-3.6.4-linux.sh
Verifying archive integrity... All good.
Uncompressing TopSpin 3.6.4 100%
NOTE: using DISPLAY=:0
Please enter root password if prompted for it.
Password:
NOTE: using DISPLAY=:0
NOTE: using TCL_LIBRARY=/tmp/selfgz1602632396/linux/tcl-8.6.9/lib/tcl8.6
NOTE: using TK_LIBRARY=/tmp/selfgz1602632396/linux/tk-8.6.9/lib/tk8.6
NOTE: running /tmp/selfgz1602632396/linux/tk-8.6.9/bin/wish8.6 -f
/tmp/selfgz1602632396/xwinstall.d/swim/lib/xwinstall.tcl --
NOTE: using log file: /tmp/install.log
```

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- Install Bruker NMR Software
- Install GPL licensed GNU GCC 4.9.1 compiler (recommended)

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- Data processing, acquisition and automation
Installs Data Processing, Acquisition and Automation Components for Spectroscopy

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- Select a spectrometer configuration to import
 - spect in /opt/topspin3.6.3

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```
*** Warning ***
```

```
Cannot import the selected spectrometer configuration (got invalid
```

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```
The IPSO computer controlling the spectrometer has to be rebooted for the
installed Diskless package to take effect.
```

```
Reboot IPSO computer automatically? [Yes]
```

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```
Error: *** Error ***
```

Installation of Spectrometer Operating System: Could not reboot the IPSO computer:
connect failed connection refused

You have to reset or reboot the IPSO computer at IP address 149.236.99.90 manually before the installed Diskless package takes effect.

To do this, either

- press the red button on the IPSO computer, or
- start TopSpin after this installation has been completed, type 'ha', open the IPSO web site, navigate to 'Administration' and click on Reboot.

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