Polígono Industrial Can Salvatella  
Avda. Salvatella, 4  
08210 Barberá del Vallés  
Barcelona - Spain  
Tel.: + 34 93 718 08 08  
Fax: + 34 93 718 23 38  
www.nirco.com

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# REPLACEMENT OF BATTERY

## Location of a nickel-metal-hydrde battery (NU-9668E/9668GC)

This unit is provided a nickel-metal-hydrde battery for the power failure warning device. The battery is located in the electrical box inside the cover on the lower left side. (Fig. 1)

The high voltage components are enclosed in the electrical box. The cover should be removed by a qualified engineer or a service personnel only to prevent the electric shock.

## Removal of nickel-metal-hydrde battery

1. Turn off the power switch and disconnect the power supply plug.
2. As shown in the Fig. 2, remove 5 screws fixing the side cover and the switch cover with a screw driver and remove the side cover.
3. Remove 4 screws fixing the electrical box cover with a screw driver. (Fig. 3)
4. Disconnect the battery connector and remove 2 screws fixing the battery mounting plate. (Fig. 4)
5. Take out the battery.
6. Follow the procedure for recycling or proper disposal.

## Handling of battery

Cover the battery terminal with an insulating tape to avoid the short circuit. Then follow the procedure for recycling or proper disposal.

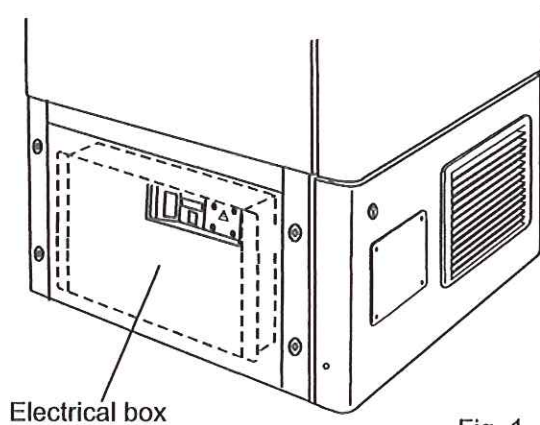
Fig. 1

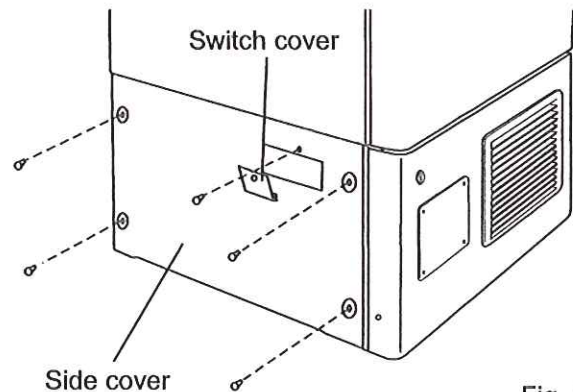
Fig. 2

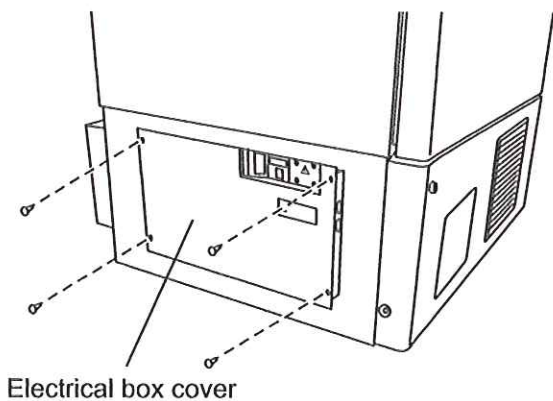
Fig. 3

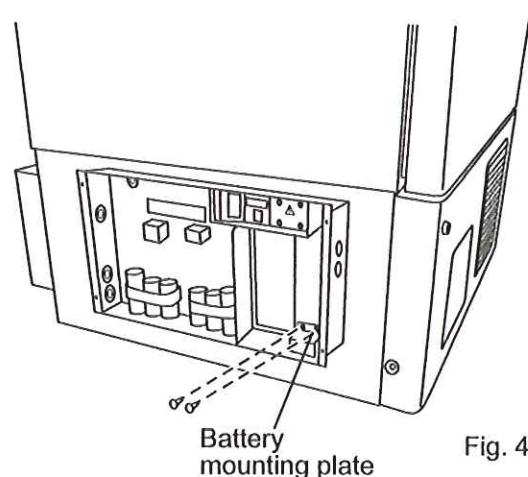
Fig. 4